

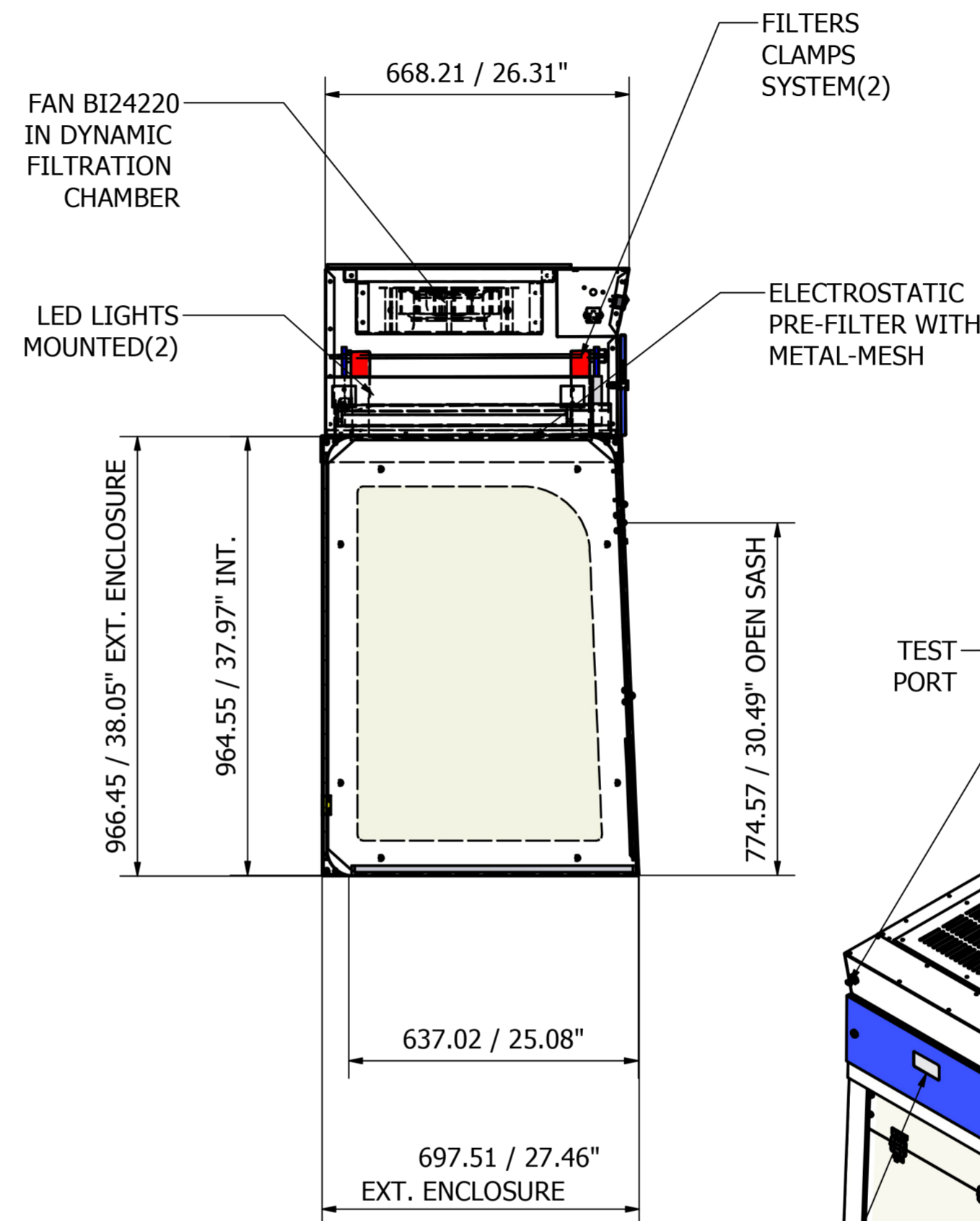


REVISION HISTORY				
ZONE	REV	DESCRIPTION	DATE	APPROVED
SIDE PANELS	1	TEMPERED GLASSES	5/18/2017	P CH
METAL HEAD	2	FAN MOD RG190	6/22/2020	P CH
METAL HEAD	3	NEW FAN MOD BI2422063TB2M-8B	6/22/2020	P CH
METAL HEAD	4	INTERNAL LED LIGHTS MOUNTED	6/22/2020	P CH
SIDE PANELS	5	CLEAR ACRYLIC ON SIDE PANELS	8/22/2022	P CH

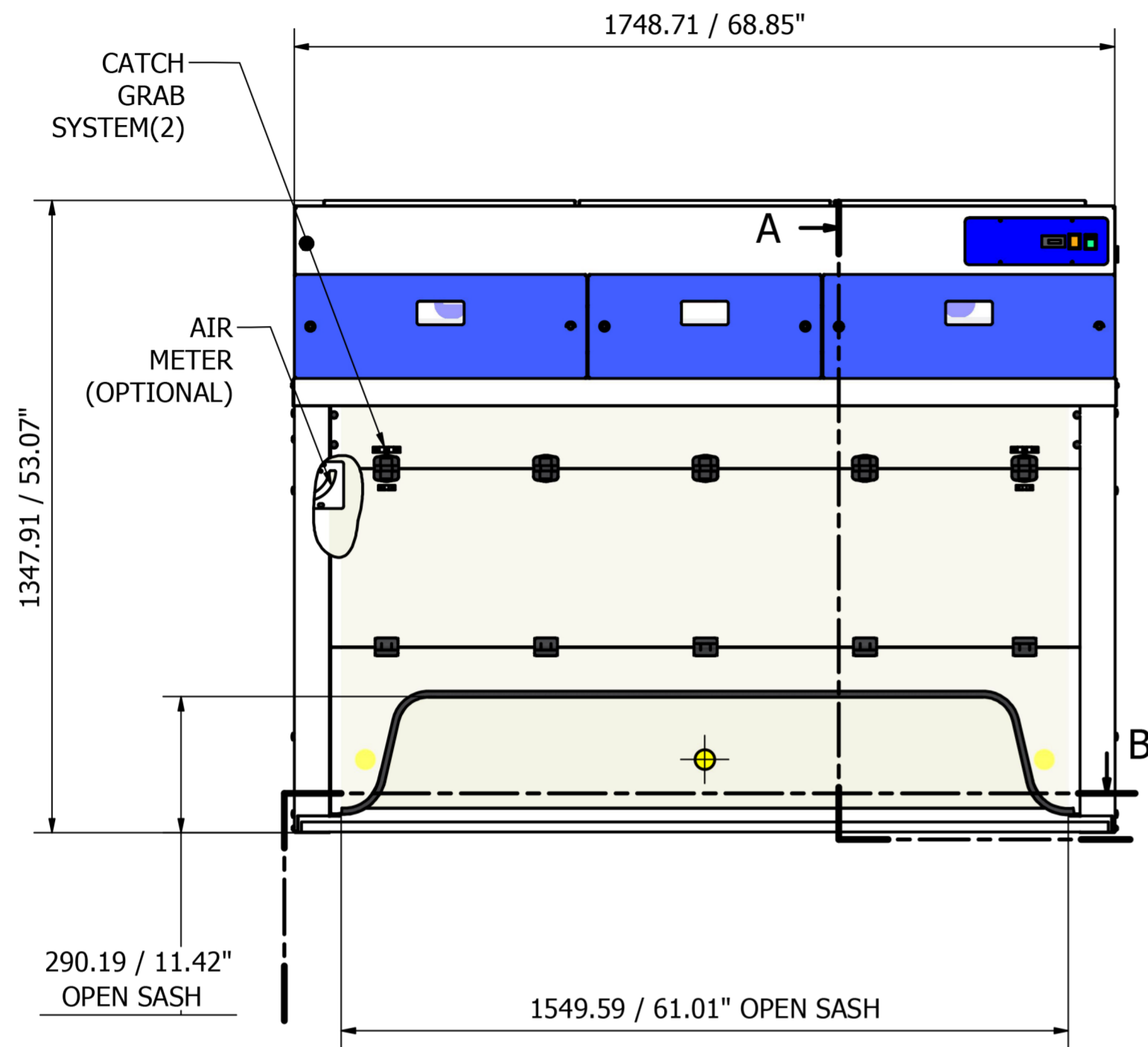

A-A ( 1 : 12 )

- NOTES:**
- STANDARD ON ALL UNITS:
  - 1.0 CONSTRUCTION: EPOXY COATED MILD STEEL SIDES PANELS AND HEAD UNIT.
  - 2.0 FRONT AND BACK PANELS MADE ON 1/4" CLEAR ACRYLIC.
  - 3.0 CLEAR ACRYLIC ON SIDE PANELS.
  - 4.0 CENTRIFUGAL FAN BI2422063TB2M-8B.
  - 5.0 POWER SUPPLY 110V 60Hz. or 220V 50Hz.
  - 6.0 MAIN SWITCH ON/OFF.
  - 7.0 MONITORING, LOW AIRFLOW.
  - 8.0 INTERNAL LED LIGHTS.
  - 9.0 W/O

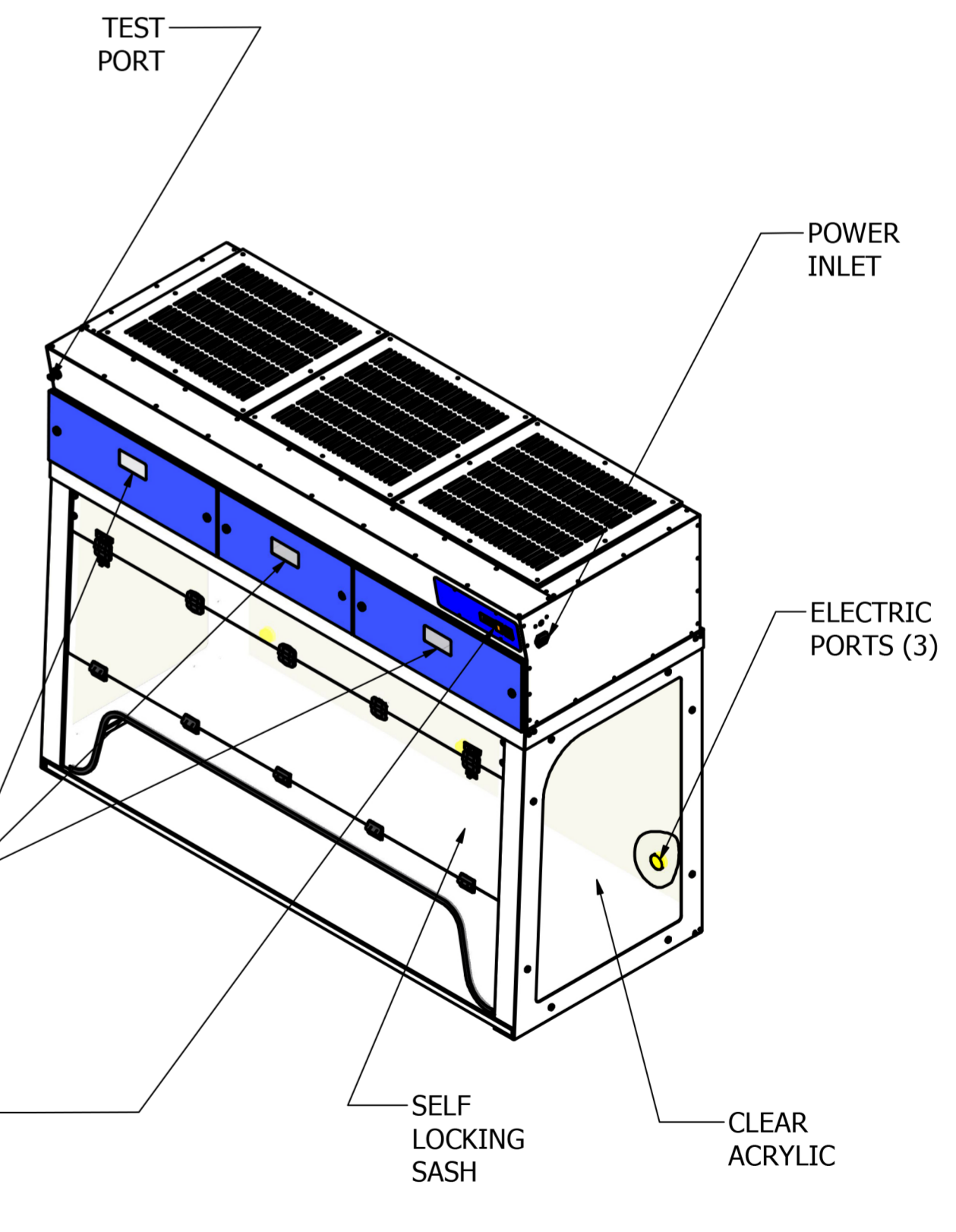

POLYPRO SPILLAGE TRAY  
25,00 x 625,00 x 1720,00  
(OPTIONAL)

B-B ( 1 : 12 )

GENERAL TOLERANCE IN MILLIMETERS	
LINEAR & ANGULAR	FORM & POSITION
ISO 2768 V	ISO 2768 L

Designed by CFG	Checked by	Approved by	Date 9/17/2012	Date
PURAIR 30XT.3F DUCTLESS FUME HOOD			Edition 5	Sheet 1 / 1
Purair 30 XT				